

In Windows10, if happen error on operating "set camera size" or "capture picture" when using coolingtech measurement software,please do as the followings:

1, After installing the software, find "CoolingTech_Microscope" by the installation file folder path ,if you choose the default path, please select the file folder "CoolingTech_Microscope" under the file folder "C:\Program Files (x86)\CoolingTech";

2, Right click on the folder "CoolingTech_Microscope" and select "properties" - > "safe" - > "Edit" button, and then click"add (D)" button, please enter "everyone"under the blank of "enter the object name to choose:" , and then click "OK" button;

3 , Please click "Everyone" under "group or user name (G):" , -->"Permission of Everyone" -->"complete control" -->"allow";

4, click on the "application" and "OK" button ,then the coolingtech measurement software can be used normally.